
Cole-Parmer®

HBC-100 & HBC-200 Series Heating Block Calibrators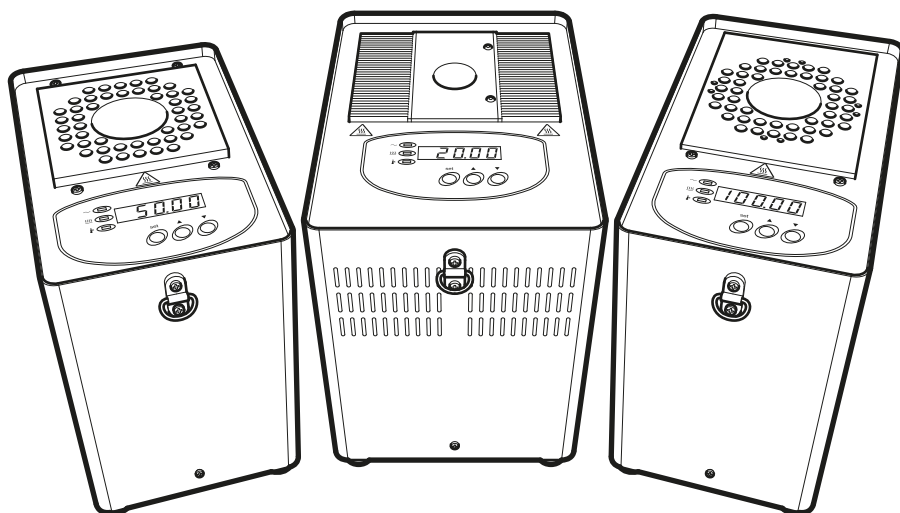

Instruction Manual
7002714-CPB Version 3

Cole-Parmer®
essentials

CONTENTS	PAGE
INTRODUCTION	3
BEFORE USE	3
SAFETY AND INSTALLATION	4
English	4
Français	6
Deutsch	8
Español	10
Italia	12
SPECIFICATION	14
OPERATION	15
Button description	15
Icon description	15
Preparation	16
Setting the operating temperature	16
Changing the display units and setting two set temperatures	16
After use	16
ADDITIONAL INFORMATION	17
User maintenance	17
Cleaning	17
Fuses	17
Fault finding	17
TECHNICAL SUPPORT AND SERVICING	17
REPLACEMENT PARTS	18
AVAILABLE INSERTS	18
SPARE PARTS	18
DECLARATION OF CONFORMITY	19

INTRODUCTION

Please read all the information in this booklet before using the unit.

The HBC-100 & HBC-200 Series Calibrators are designed to provide safe and convenient calibration of a wide range of thermal sensors. It features fast heat up times, with accuracy and repeatability.

Applications

The HBC unit can calibrate temperature probes without the need to return them to a specialist laboratory. To ensure accuracy the unit must be operated in an environment with a stable ambient temperature.

The thermal sensors are placed in a well in the temperature controlled block. A number of inserts with different dimension wells are available to match standard probe sizes: these are detailed towards the back of this manual.

Maintaining a set temperature

The required temperature is set on the calibrator and the operation of the probe is checked. ukAS certificate can be provided, contact your dealer for details.

BEFORE USE

Before using the Calibrator please make sure you have read this manual carefully. If there is any doubt relating to the proper use of this equipment, the staff at Antylia Scientific Ltd. or your supplier will be happy to assist you.

Unpacking

When unpacking the unit please ensure that the following have been removed from the packaging:

HBC-100 or HBC-200 Series Block Calibrator

Extraction tool

Mains cables (UK and EU or US)

Operator's manual

The user is advised to keep the original packaging in case the instrument ever needs to be returned for service or repair. Antylia Scientific Ltd. accepts no responsibility for damage incurred unless the unit is correctly packed and transported in its original packaging.

SAFETY AND INSTALLATION

Please read all the information in this manual before using the Calibrator.

WARNING

HIGH TEMPERATURES ARE DANGEROUS: they can cause serious burns to operators and ignite combustible material.

Cole-Parmer have taken great care in the design of these units to protect operators from hazards, but operators should pay attention to the following points:

- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS.
- DO NOT use combustible substances near hot objects.
- DO NOT operate the instrument in the vicinity of inflammable liquids or gases.
- DO NOT place any liquid directly into the instrument.

Operator Safety

All operators of Cole-Parmer equipment must have available the relevant literature needed to ensure their safety. It is important that only suitably trained personnel operate this equipment, in accordance with the instructions contained in this manual and with general safety standards and procedures. If the equipment is used in a manner not specified by Antylia Scientific Ltd. the protection provided by the equipment to the operator may be impaired.

All Cole-Parmer instruments are designed to conform to international safety requirements and are fitted with an over-temperature cut-out.

If a safety problem should be encountered, switch off the unit at the mains socket and remove the plug from the electricity supply.

Installation

The unit should be carried using both hands. Never move or carry the instrument when in use or connected to the mains electricity supply.

1. All Cole-Parmer instruments are supplied with a power cable; this may be integral or plug-in.

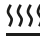
2.  Before connecting the instrument to the mains electricity supply, check the voltage against the rating plate (located on the back of the unit). **Please note that the unit must be earthed to ensure proper electrical safety.** Connect the mains cable to a suitable plug according to the table below.

Connections	220V-240V, 50/60Hz	110V-120V, 50/60Hz
Live	Brown	Black
Neutral	Blue	White
Earth	Green/yellow	Green

Note that units marked 230V on the rating plate work at 220V; units marked 120V work at 110V. In both cases, however, the heating rate will degrade by approximately 8%.

3. Place the unit on a suitable flat bench or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction.

4. Plug the mains cable into the socket on the back of the instrument.
5. Note that the following symbols may be next to the indicator lamps on the front panel of the units and have the following meanings:

-  the power indicator
-  the heater indicator
-  the over-temperature indicator

6. Switch on the instrument. Symbols on or near the power switch of the unit have the following meanings:
 - I Mains Switch On
 - O Mains Switch Off

Replacement cable

Should the mains lead need replacement, a cable of 1mm² of harmonized code H05VV-F connected to an IEC320 plug should be used. **IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN.**

Environmental Conditions

The Calibrator is designed to operate under the following conditions:

- Indoor use
- Ambient temperature range +5°C to +40°C
- Altitude to 2000m
- Relative humidity up to 95%
- Mains supply fluctuations not exceeding 10%
- Over voltage category II IEC 60364-4-443
- Pollution degree 2
- Use with a minimum distance all-around of 200mm from walls or other items

Note: The control specifications are quoted at an ambient temperature of 20°C. The specification may deteriorate outside an ambient temperature of between 10°C and 30°C.

The unit has been tested for radio frequency interference and is certified under EN61326.

Guarantee

The unit is guaranteed against any defect in material or workmanship for the period specified on the enclosed guarantee card. This period is effective from the date of purchase; within this period all defective parts will be replaced free of charge provided that the defect is not the result of an accident, misuse or negligence. Servicing under this guarantee should be obtained from the supplier of the instrument.

This manual has been prepared for the convenience of Cole-Parmer customers and nothing in this manual shall be taken as a warranty, condition or representation concerning the description, merchantability, fitness for purpose or otherwise of the unit or components.

Notwithstanding the description and specification(s) of the instruments contained in the operator's manual, Cole-Parmer reserves the right to make such changes as it sees fit to the instruments or to any of the components.

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SPECIFICATION


HBC-100-140F Specification

The figures stated in the specification are as measured 0-50 mm from the base of the well of the insert, using an SPRT, at the time of calibration.

Minimum temperature	*45°C below ambient Typically -20°C in an ambient of 25°C
Maximum temperature	140°C
Temperature accuracy	±0.3°C
Temperature uniformity	±0.2°C
Temperature stability after 10 minutes	±0.05°C
Display resolution	0.01°C
Set point display resolution	0.1°C
Heat up rate 5 minutes	20°C to 100°C
Cool down rate 9 minutes	100°C to 0°C
Fan cooling Automatic	Automatic
Dimensions Height x Width x Depth	273mm x 207mm x 289mm
Large well in the block:	Depth / Diameter 130mm / 38.2mm
Weight	11Kg

Options

- Central Well for inserts, see the list of inserts
- Full bi-directional RS232 Comms Port 9way D type plus software and cable
- Soft carry case
- Hard carry case
- UKAS calibration certificate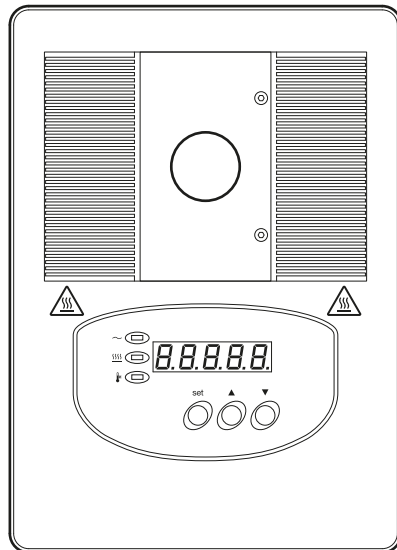

*Although temperature can be set to -40°C this can only be achieved in an ambient of 5°C.

HBC-200-425F Specification

The figures stated in the specification are as measured 0-50 mm from the base of the well of the insert, using an SPRT, at the time of calibration.

Minimum temperature	20°C above ambient
Maximum temperature	425°C
Temperature accuracy	±0.3°C
Temperature uniformity	±0.2°C
Temperature stability at 200 after 10 minutes	±0.03°C
Temperature stability at 425 after 10 minutes	±0.05°C
Display resolution	0.01°C
Set point display resolution	0.1°C
Heat up 12 minutes	20°C to 400°C
Cool down 21 minutes	400°C to 100°C
Fan cooling Automatic	Automatic
Dimensions Height x Width x Depth	270mm x 170mm x 255mm
Large well in the block:	Depth / Diameter 130mm / 38.2mm
Weight	6.3Kg

Options


Central Well for inserts, see the list of inserts

Full bi-directional RS232 Comms Port 9way D type plus software and cable

Soft carry case

Hard carry case

UKAS calibration certificate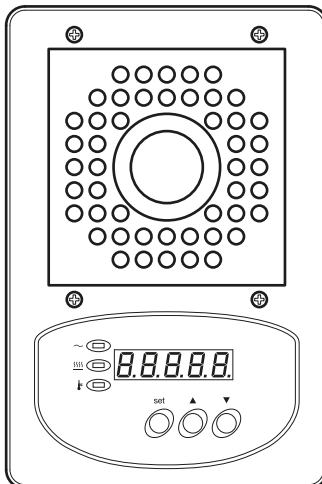

HBC-200-650F Specification

The figures stated in the specification are as measured 0-50 mm from the base of the well of the insert, using an SPRT, at the time of calibration.

Minimum temperature	25°C above ambient
Maximum temperature	650°C
Temperature accuracy	±0.4°C
Temperature uniformity	±1°C
Temperature stability after 10 minutes	±0.09°C
Display resolution	0.01°C
Set point display resolution	0.1°C
Heat up 35 minutes	20°C to 600°C
Cool down 30 minutes	600°C to 200°C
Fan cooling Automatic	Automatic
Dimensions Height x Width x Depth	278mm x 170mm x 300mm
Large well in the block:	Depth / Diameter 152mm / 38.2mm
Weight	9.6Kg

Options


Central Well for inserts, see the list of inserts

Full bi-directional RS232 Comms Port 9way D type plus software and cable

Soft carry case

Hard carry case


UKAS calibration certificate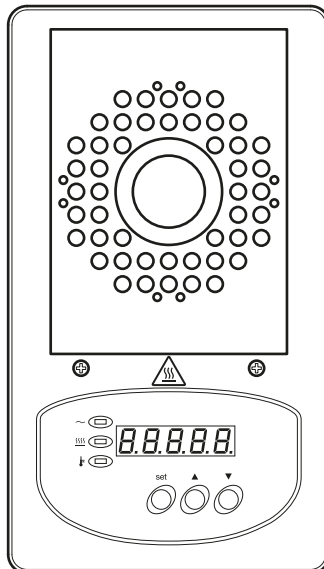


OPERATION

When You Switch On

When you first switch on, the display will show the edition of the software which your unit uses. For example software issue "3.13" would be shown as follows: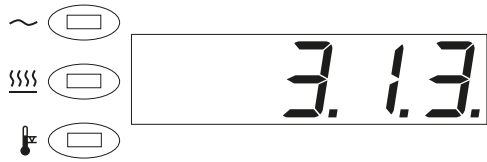

It will display this for 1 second, then the actual temperature of the block will be indicated.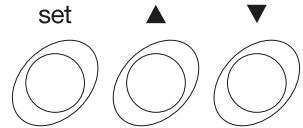


The Front Panel Controls

The front panel controls consist of three buttons for controlling the display, a five digit LED display and three indicators.


The SET temperature Button

The SET temperature button displays the set temperature when pressed.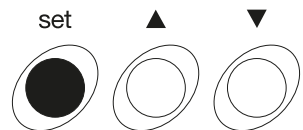

The UP ARROW Button

When the SET temperature button is held down and the UP ARROW button is pressed, the set temperature is increased.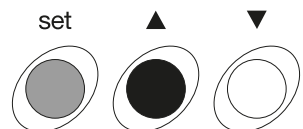


The DOWN ARROW Button

When the SET temperature button is held down and the DOWN ARROW button is pressed, the set temperature is decreased.

Speed of Change of Set Temperature

Each press of the UP ARROW or DOWN ARROW buttons will increase or decrease the set temperature by 0.1°C. If the buttons are held down the temperature change will accelerate to 5° per second.

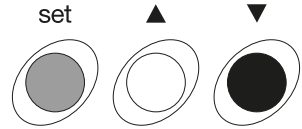Power Indicator

The top indicator shows that there is power to the unit

Power to the Block Indicator

The next indicator shows when there is power to the block.

When the set temperature is different to the block temperature this indicator will light. If the light is on continuously the block is getting constant power. The only exception is described under Over-Temperature Indicator. As the temperature approaches the set temperature the indicator will flash. When set temperature is reached the indicator will stay on for shorter periods.

Sensor fault Indicator

If there should, for any reason, be a sensor fault, the bottom indicator will light. The power to the block will have been switched off and the unit will begin to return to ambient even if the power light is on (the light staying on or not depends on which circuit has sensed a fault).

Setting the units for temperature

The unit will normally be set to display in °C, if you require it to work and display in °F press both the UP ARROW and DOWN ARROW buttons at the same time. To convert back, press both buttons again.

Setting the Operating Temperature

1. To display the set temperature on the digital display, press and hold the SET Temperature button. To adjust the set temperature, press the SET Temperature button and hold it while pressing the UP or DOWN buttons. When the SET Temperature button is released, the measured temperature is displayed (in degrees Celsius).
2. Power is switched on to the block (and the indicator comes on) if the set temperature is higher or lower than the block temperature.
3. When the measured temperature approaches the set temperature, the "power to the block" indicator will begin to flash. As the measured temperature stabilises the indicator will stay on for shorter periods.
4. There will be a time-lag between the block and the insert achieving the set temperature due to thermal contact between them. The display may reach temperature in 2 or 3 minutes while the insert may take 5 minutes or more to reach temperature.

Overtemperature cut-out

The unit is fitted with an overtemperature cut-out. It will operate if a fault develops and the unit exceeds its maximum operating temperature. The cutout will remove power to the heaters or peltiers and return the unit to ambient temperature.

If the cut-out trips, the unit needs to be turned off and then turned on again to reset it.

If the unit is switched off and then switched on again quickly this may cause the overtemperature cut-out to trip. If this occurs, turn the unit off, wait 5 seconds or more then switch it on again.

Repeated tripping of the cut-out may indicate a fault and you should contact your supplier for more information.

After Use

When you have finished heating samples, remember that parts of the unit inserts and associated accessories may be very hot. Take the precautions listed earlier. We recommend that the inserts should be allowed to cool to 70°C before being removed from the HBC unit. They will still have to be handled with care


If you need to remove an insert while it is hot, fit the extractor tool into the locating holes and lift the insert out carefully. Never leave the extractor tool in the insert while it is being used in the HBC unit.

RS232 Serial Interface (where applicable)

The calibrator may send data logging information to a PC or compatible computer by connecting the unit and the PC via an RS232 cable, and installing the "Calsoft" software, available free of charge from our web-site.

Ensure that there is a connection made before the calibrator is switched on.

The RS232 cable must be fitted to both the unit and the PC before either unit is powered up, otherwise, data integrity cannot be guaranteed. Once the cable is fitted, it does not matter which unit is powered up first.

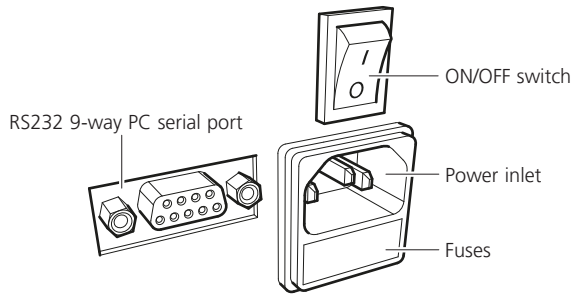The following tables indicate the cable specifications for a 9-way PC serial port:

Calibrator		PC
9-way female D type		9-way female D type
pin	signal	pin
Case	F.GND	Case
3	TxD	3
2	RxD	2
7	RTS	7
8	CTS	8
6	DSR	6
1	CD	1
4	DTR	4
5	S.GND	5

Sending data to the PC

The procedure for sending data to the PC is described in the instructions supplied with the software. The PC must be properly connected by the RS232 cable and running Calsoft to accept or display data.

ADDITIONAL INFORMATION

User Maintenance

NOTE THAT THIS EQUIPMENT SHOULD ONLY BE DISMANTLED BY PROPERLY TRAINED PERSONNEL. REMOVING THE FRONT OR REAR PANELS EXPOSES POTENTIALLY LETHAL MAINS VOLTAGES. THERE ARE NO OPERATOR MAINTAINABLE PARTS WITHIN THE EQUIPMENT.

In the unlikely event that you experience any problems with your unit which cannot easily be remedied, you should contact your supplier and return the unit if necessary. Please include any details of the fault observed and remember to return the unit in its original packing. Cole-Parmer will accept no responsibility for any damage to units that are improperly packed for shipment. If in doubt, contact your supplier. See the Decontamination Certificate supplied with your unit.

Cleaning

Before cleaning your unit, ALWAYS disconnect it from the power supply and allow it to cool below 50° C. Your unit can be cleaned with a cloth dipped in water or ethanol (methanol can also be used). No part of the case or cover should be immersed in the solvents. Do not use acetone or abrasive cleaners.

Before using any cleaning or decontamination method except those recommended in this manual, the responsible body should check with Cole-Parmer that the proposed method will not damage the equipment.

Fuses

Your unit is protected by two fuses. They should only be changed by suitably qualified personnel. If the fuses blow persistently, a serious fault is indicated and you may need to return the unit to your supplier for repair.

Fault Finding

HELP display: If HELP is displayed there is a problem with the internal sensing device. The calibrator should be returned to your supplier for repair.

Over-temperature cut-out: If the over-temperature cut-out trips it can be reset by switching the unit off, waiting a few seconds, then swiftly it on again. If it repeatedly trips, please contact your local supplier.

Should you have any problems with your Calibrator which cannot be easily remedied, you should contact your supplier and return the unit if necessary. Please include details of the fault observed and remember to return the unit in its original packing. Antylia Scientific Ltd. accepts no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier, giving the full serial number of the unit.

TECHNICAL SUPPORT AND SERVICING

If you require further technical or application assistance please contact:

Email: cptechsupport@antylia.com

Phone: +44 (0)1785 810433

For servicing information please contact:

Antylia Scientific Ltd. Service Department, Beacon Road, Stone, Staffordshire, ST15 0SA,

Email: cpsservice@antylia.com.

Phone: +44 (0)1785 810475

We are continually striving to improve our products and software. If you have any comments and suggestions on how we can do things better please send them to us at: cptechsupport@antylia.com

REPLACEMENT PARTS

Part No.	Description
6101058	Insert extractor
FCAB10UK	Mains cable, 230V UK units
FCABLEEU	Mains cable, 230V units
FCABLEUS	Mains cable, 120/100V units

ACCESSORIES

Part No.	Description
FCAL232	CALSOFT PC software kit (includes RS232 cable) Supplied with models with RS232 connection
FCALFCSE	Soft carrying case
FCALFCSH	Hard Carrying case
FDB00CP	Cooling probe 425 & 650 requires an insert with a 10mm diameter or larger hole
6103890	RS232 Cable only

INSERTS

Inserts are made from aluminium or aluminium-bronze and must be ordered separately from the calibrator. Each insert is stamped for identification.

		HBC-100-140F and HBC-200-425F Inserts				HBC-200-650F Inserts	
		Aluminium Inserts: Immersion depth 114.3 mm or 4.5"				Aluminium Bronze Inserts: Immersion depth 152.4 mm or 6"	
		Part No.	Diameter			Part No.	Diameter
Metric	FINSALA	5 off 6mm		Metric	FINSABA	5 off 6mm	
	FINSALB	1 off each 10mm, 8mm, 6mm, 4.5mm, 3mm			FINSABB	1 off each 10mm, 8mm, 6mm, 4.5mm, 3mm	
	FINSALC	2 off 6mm, 2 off 10mm			FINSABC	2 off 6mm, 2 off 10mm	
	FINSALD	2 off 6mm, 2 off 12mm			FINSABD	2 off 6mm, 2 off 12mm	
	FINSALE	1 off 6mm			FINSABE	1 off 6mm	
	FINSALZ	1 off 20mm			FINSABZ	1 off 20mm	
Imperial	FINSALF	5 off 1/4"		Imperial	FINSABF	5 off 1/4"	
	FINSALG	1 off each 3/8", 5/16", 1/4", 3/16", 1/8"			FINSABG	1 off each 3/8", 5/16", 1/4", 3/16", 1/8"	
	FINSALH	2 off 1/4", 2 off 3/8"			FINSABH	2 off 1/4", 2 off 3/8"	
	FINSALI	2 off 1/4", 2 off 1/2"			FINSABI	2 off 1/4", 2 off 1/2"	
	FINSALJ	1 off 1/4"			FINSABJ	1 off 1/4"	
	FINSALL	1 off 9/16"			FINSABL	1 off 9/16"	
	FINSALM	1 off 5/8"			FINSABM	1 off 5/8"	
	FINSALN	1 off 3/4"			FINSABN	1 off 3/4"	
FINSALO	1 off 11/16"		FINSABO	1 off 11/16"			
Blank	FINSALK	BLANK (A blank insert may be drilled to suit your application)		Blank	FINSABK	BLANK (A blank insert may be drilled to suit your application)	

This product meets the applicable CE Directives and UKCA Legislation for radio frequency interference and may be expected not to interfere with, or be affected by, other equipment with similar qualifications. We cannot be sure that other equipment used in its vicinity will meet these standards and so we cannot guarantee

that interference will not occur in practise. Where there is a possibility that injury, damage or loss might occur if equipment malfunctions due to radio frequency interference, or for general advise before use, contact the manufacturer.

Declaration of Conformity is available to view online at www.coleparmer.com

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Cole-Parmer®
essentials

Antylia Scientific Ltd.

Beacon Road,
Stone,
Staffordshire,
ST15 0SA,
United Kingdom